

Whatman™ Nuclepore™ Track-Etched Membranes

Nuclepore Polycarbonate
Nuclepore Polyester

Product Information sheet

Warning

For research use only.

Not recommended or intended for diagnosis of disease in humans or animals.

Do not use internally or externally in humans or animals.

Membrane Information Guide

	Nuclepore (PC) Polycarbonate	Nuclepore (PE) Polyester
Orientation for Use	Use shiny side up* Packaged shiny side down.	Use as packaged. Both sides are the same.
Hydrophilic	Yes	Yes
Wetting Agent	PVP (Polyvinylpyrrolidone) or PVP free	PVP (Polyvinylpyrrolidone) or PVP free
Autoclaving	Autoclavable 121°C	Autoclavable 121°C
Retractive Index	1.584, 1.625	1.55-1.66
Solvents for Dissolving Membranes	N-methyl-2-pyrrolidone Methylene chloride Chloroform	Hexafluoroisopropanol M-Cresol O-Chlorophenol Phenol tetrachloroethane (1:1)
Chemical Compatibility	See catalog Chemical Compatibility Chart	See catalog Chemical Compatibility Chart

*For microscopy.

Handling Tips:

1. To avoid contamination, use forceps when handling the membrane.
2. Do not touch the membrane with your fingers.
3. Remove the separator paper before using the membrane.

Limited Warranty Disclaimer

OUR MEMBRANE FILTRATION PRODUCTS ARE SOLD FOR INDUSTRIAL AND GENERAL LABORATORY USE ONLY. NOT INTENDED AND NOT TESTED FOR MEDICAL OR THERAPEUTIC USE. We warrant to purchasers that for one year from the date of purchase we will replace any of our membrane filtration products which are defective in materials and workmanship. This is the only warranty and ANY OTHER WARRANTIES (INCLUDING ANY IMPLIED WARRANTIES OF MERCHANTABILITY AND FITNESS FOR A PARTICULAR PURPOSE) ARE EXCLUDED. The Warranty does not apply to membrane filtration products which are altered, broken or damaged by misuse or failure to follow the provided instructions.

Legal.

GE, imagination at work and GE Monogram are trademarks of General Electric Company.

Whatman and Nuclepore are trademarks of GE Healthcare companies.

© 2007–2013 General Electric Company – All rights reserved.
Previously published Dec 2007.

All goods and services are sold subject to the terms and conditions of sale of the company within GE Healthcare which supplies them. A copy of these terms and conditions is available on request. Contact your local GE Healthcare representative for the most current information.

For your local office contact information, visit
www.gelifesciences.com/contact

GE Healthcare UK Limited
Amersham Place
Little Chalfont, Buckinghamshire,
HP7 9NA, UK

<http://www.gelifesciences.com>

GE Healthcare offices:

GE Healthcare Bio-Sciences AB
Björkgatan 30, 751 84 Uppsala,
Sweden

GE Healthcare Europe GmbH
Munzinger Strasse 5, D-79111 Freiburg,
Germany

GE Healthcare Bio-Sciences Corp.
800 Centennial Avenue, P.O. Box 1327,
Piscataway, NJ 08855-1327,
USA

GE Healthcare Japan Corporation
Sanken Bldg. 3-25-1, Hyakunincho,
Shinjuku-ku, Tokyo 169-0073,
Japan



imagination at work